

(F) Microscope de poche X30 avec mini lampe

Attention : N'observez pas le soleil ou une forte source de lumière avec le microscope ce qui pourrait nuire sérieusement à la vue. Ne convient pas aux enfants de moins de huit ans.

Caractéristiques techniques :


- Grossissement x 30
- Ampoule 2.2v 250 mA
- Fonctionne avec deux piles de 1,5 volts (non incluses) type AAA
- Corps en plastique abs
- Lentilles en verre

Utilisation :


- Positionner le microscope verticalement dans le support en plexiglas.
- Positionner la lampe en biais de l'autre côté du support.
- Allumer la lampe avec l'interrupteur situé sur le côté de celle-ci.
- Regardez dans le microscope et régler l'image à l'aide du bouton 'FOCUS' du microscope. On peut aussi l'utiliser sans lampe et support.

Mise en place des piles :

- Appuyez sur les deux clips situés sur chaque face de la lampe.
- Retirer le capuchon gris.
- Placer les piles en respectant la polarité indiquée sur la lampe.
- Replacer le capuchon gris.
- Vérifier le fonctionnement avec l'interrupteur.

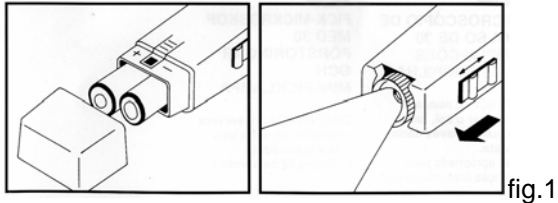

Remplacement de la lampe :

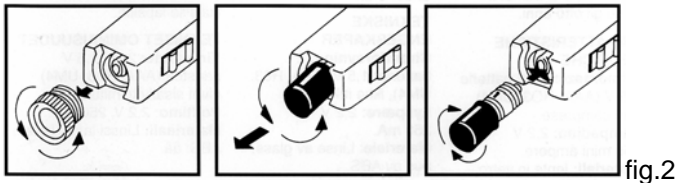

- Devisez la lentille.
- Utilisez l'embout noir pour deviser la lampe en faisant coïncider les deux ergots.
- Replacer la nouvelle lampe de même façon.

(GB) Pocket Microscope X30 with mini lamp

Caution: Do not observe the sun or a strong source of light with the microscope what could harm the sight seriously. Not suitable for children under 8 years.

Design features:

- Magnification X 30
- Bulb 2.2v 250 mA
- Function with two batteries 1,5 volts (not included) standard AAA
- Plastic ABS body
- Lenses out of glass

Use:

- Put in vertically position microscope in the plexiglass support.
- Put the lamp close to the microscope in the support in 45° position.
- Light the lamp with the switch located on the body.
- Look in the microscope and to regulate the image with ' FOCUS' button of the microscope
- It can be use without support and lamp.

Installation of batteries (fig. 1):

- Press on the two clips located on each face of the lamp.
- Pull up the grey cap.
- Place the batteries by respecting the polarity indicated on the lamp.
- Replace the grey cap.
- Check operation with the switch.

Replacement of the bulb (fig. 2):

- Unscrew the lens.
- Use the black end for unscrews the bulb.
- Replace the new bulb in the same way.